

The Seedburo Sheller provides a rapid, easy and efficient method of shelling. Utilizing only one piece of equipment it will accommodate both wet and dry rice samples for all varieties of rough rice. The Seedburo Sheller is designed to comply with service requirements of the Federal/State Rice Inspection Service. Quality construction assures years of service. Safety features prevent damage by small rocks or metal objects, which accidentally enter the hopper. The large hopper holds over 2 lbs. of rough rice. This unit will shell 1 lb. per minute in continuous operation. The dust collector is used to save hulls from each sample for further analysis.



**Sheller:** Net Wt. 124 lbs., Ship Wt. 166 lbs, Ship Dims. 33" L x 21" W x 31" H  
**Dust Collector:** Net Wt. 16 lbs., Ship Wt. 35 lbs, Ship Dims. 16" L x 15" W x 28" H

Sample Sheller with Dust Collector, 1/3 HP, 115V, 60Hz .....	<b>No. 64-115-60-WDC</b>
Sample Sheller with Dust Collector, 1/3 HP, 220V, 50Hz .....	<b>No. 64-220-50-WDC</b>
Sample Sheller without Dust Collector, 1/3 HP, 115V, 60Hz .....	<b>No. 64-115-60-WODC</b>
Sample Sheller without Dust Collector, 1/3 HP, 220V, 50Hz .....	<b>No. 64-220-50-WODC</b>
Crating for Sheller .....	<b>No. 64/DOM</b>

## RICE SIZING DEVICE

### Here's How It Works:

First, insert two plates of the same size into the machine, then feed a standard sample of 1000 grams of cleaned shelled, milled rice onto the top screen. When the entire sample has passed over the plates, remove and empty the rice that remains on the plates into a separate container. Change to the next set of plates and repeat the above steps, until all necessary separations have been made.

Because the total sample was 1000 grams, the percentage of each portion is obtained by weighing it and simply moving the decimal point one place. The percentage of whole rice is taken from the weight of the rice which remains after the last set of plates have been used.

**Rice Sizing Device:** Net Wt. 110 lbs, Ship Wt. 160 lbs. Ship Wt with Set of 10 Plates 200 lbs, Ship Dims. 36" L x 35" W x 20" H  
**Single Plate:** Net wt.3 lbs., Ship Wt. 5 lbs, Ship Dims 25" L x 11" W x 4" H.

Rice Device, less plates, 115V, 60Hz .....	<b>No. 61-115-60</b>
Rice Device, less plates, 230V, 60Hz .....	<b>No. 61-230-60</b>
Rice Device, less plates, 110V, 50Hz .....	<b>No. 61-110-50</b>
Rice Device, less plates, 220V, 50Hz .....	<b>No. 61-220-50</b>
Crating, Wooden Box .....	<b>No. 61/DOM</b>

### Standard set of 10 plates consists of:

Plates perforated with round holes, centers staggered. Two of each size: 5/64" (centers 5/32") 6/64" (centers 5/32"), .135" (centers 3/16"), 10/64" (centers ¼"), 12/64" (centers ¼").

Standard set of 10 plates .....	<b>No. 61SET</b>
12/64" Plate .....	<b>No. 61-12</b>
10/64" Plate .....	<b>No. 61-10</b>
6/64" Plate .....	<b>No. 61-6</b>
5/64" Plate .....	<b>No. 61-5</b>
.135" Plate .....	<b>No. 61-135</b>
13/64" Plate .....	<b>No. 61-13</b>

